### **Product Features**

Low-noise, battery-based design with  $815pA/\sqrt{Hz}$  noise density

10ppm stability over 30 minutes

Constant current and constant power operating modes

Dual current ranges: 200mA and 500mA floating output

1MHz external modulation bandwidth

Programmable internal ramp generator

Built-in battery charger

The LDX-3620 is a battery-powered, ultra-low noise current source, optimized for narrow linewidth or stable wavelength laser diode applications. The instrument provides a dual-current range with a 4.5-digit display. It offers high stability performance, constant current or power operation, internal or external modulation capability, and comes with a built-in battery charger.



Ultra-Low Noise Current Source



High Stability Performance with Flexible Modulation Capability



# LDX 3620

# www.yawawetigeni Current Source

### Ultra Low Noise and High Stability

Careful component selection and circuit board

CS\_location liver unprecedented low noise levels.

Current noise density is only 815pA/√Hz.

Wideband current noise is only 850nA rms. The unique design also delivers stable performance that outperforms conventional current sources.

Even over periods of tens of minutes, output is constant to within 10ppm. Such attention to noise and drift figures is critical for applications such as laboratory work in coherent communications or atomic spectroscopy.

### **Internal or External Modulation**

The LDX-3620 offers two ways to modulate the output. One, the built-in programmable ramp generator lets you make an L/I curve or dither the wavelength without introducing noise from an external signal generator. In addition, a trigger signal on the rear of the instrument permits oscilloscope triggering at ramp start-up. Two, an external signal of up to 1MHz-bandwidth, AC- or DC-coupled, and input through the front panel BNC connector modulates the output current. In all modes of operation, the modulation signal is summed with the constant current output.

### Laser Diode Protection


The instrument's slow-start circuit ensures a gentle laser turn-on. For maximum protection, this circuit also keeps the output shorted during power-up and when changing lasers. The LDX-3620 also employs a redundant current limit that may be set using the recessed front panel knob. This safety feature provides extremely fast and effective current limiting under all modes of operation.

### **Built-In Battery with Charger**

The LDX-3620's internal batteries provide the benefits of low noise without the normal drawbacks of battery-powered instruments. The instrument's lead-acid batteries offer high charge retention and long operating times. A battery charger is built into the instrument, permitting instrument operation even while the batteries are being charged.

### **Low Battery Protection**

The 3620 has two low battery circuits. The first simply displays "LO BAT" when the batteries need recharging. The second shuts the instrument down safely if the battery charge gets dangerously low.

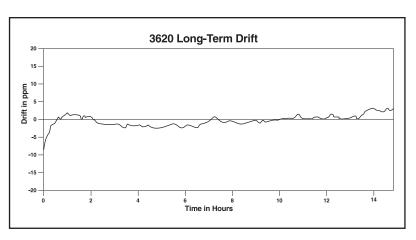

The 3620 performs with -10ppm drift (after 30 minutes) and 0.2% constant power stability.

# LDX 3620

Ultra-Low Noise Current Source



# Ultra-low noise current source that's optimized for your laser diode application.

### **Constant Current Operation**

When operated in constant current mode, all unused circuitry is switched out, allowing measured output noise figures to approach the theoretical limits for the control circuit. The 4.5-digit liquid crystal display, when used in conjunction with the coarse and fine output controls, allows precise control of the current to your laser.

### **Constant Power Operation**

When in constant power mode, the photocurrent from a back facet monitor photodiode or a front facet power monitor may be used to maintain constant output power. Controls on the rear of the instrument allow the gain to be changed depending on the magnitude of the photocurrent.

# www.yayawetiganics.(dayayayidz bandwidth) **Current Source**

### Specifications<sup>1</sup>

OUTPUT	0-200 mA	0-500mA
Output Current:2	0-200mA	0-500mA
Compliance Voltage:	<u>&gt;</u> 5V	<u>≥</u> 4V
Noise and Ripple		
(54 1-4-400 A) 1- 1- 1- 1- 1- 1- 1- 1- 1- 1- 1- 1- 1-		

<2μA rms ≤850nA rms AC line operation: <3µA rms <8µA rms (5Hz-10kHz bandwidth) ≤500nA rms Battery operation: ≤100nA rms AC line operation: <6.5µA rms <2µA rms Noise Density  $5.2 \text{nA}/\sqrt{\text{Hz}}$ 11.5nA/ $\sqrt{\text{Hz}}$ 50/60Hz:

Battery Operation: 815pA/ $\sqrt{\text{Hz}}$  $2.6 \text{nA}/\sqrt{\text{Hz}}$ 1kHz: 25kHz: 315pA/√Hz 795pA/√Hz

Stability 10-20 seconds: <0.1ppm <0.1ppm 3-5 minutes: <1ppm <1ppm 10-30 minutes: <10ppm <10ppm ≤10ppm/°C Temperature Coefficient: <10ppm/°C Transients:  $\leq 10 \mu A$  $\leq 10 \mu A$ Shielded 9-pin Shielded 9-pin Output Connector:

D-sub on panel D-sub on panel

**CURRENT LIMIT** 

10-500mA 10-500mA Range: Accuracy: ±5mA ±5mA

DISPLAY

Connectors:

4.5-digit LCD 4.5-digit LCD Type: Range 0-199.99 0-500.0 Resolution: 0.01mA 0.1mA ±0.05mA ±0.6mA Accuracy: LASER diode LASER diode Reads: current in mA current in mA **MONITOR MONITOR** photodiode photodiode current in mA current in mA

PHOTODIODE FEEDBACK

Current input Current input Input Type: from external from external photodiode photodiode Range: 20µA to 2mA 20µA to 2mA for full scale for full scale

output output Rear panel

Rear panel isolated BNC isolated BNC iack or 9-pin jack or 9-pin rear panel rear panel D-sub D-sub connector connector

INTERNAL RAMP GENERATOR

Period: Adjustable Adjustable 50ms to 100s 50ms to 100s

Span Adjustment: 0-200/ 0-200/ 0-500mA 0-500mA



P.O. Box 6310, Bozeman, MT 59771 • FAX: 406-586-9405

www.ilxlightwave.com

Ramp Trigger: Ramp start Ramp start trigger output trigger output

> for oscilloscope; for oscilloscope; optically optically isolated open isolated open collector TTL collector TTL

output output

Ramp Flyback Time: 700µs, 700µs, approximately approximately

EXTERNAL MODULATION INPUT

Bandwidth (3dB point)

AC Coupled: 100Hz to 1MHz 100Hz to 1MHz DC Coupled: DC to 1MHz DC to 1MHz 250mA/V Transfer Function: 100mA/V Connector: Isolated front Isolated front panel BNC panel BNC

**GENERAL** 

AC Power Input Voltage

> Range: 100-125 or 100-125 or 210-250VAC 210-250VAC

Line Frequency: 50/60Hz 50/60Hz

Battery

+12 volt supply: 4.5A-hour sealed 4.5A-hour sealed

lead-acid battery lead-acid battery (standard) (standard)

-12 volt supply: 1.2A-hour sealed 1.2A-hour sealed lead-acid battery lead-acid battery

Temperature Range

0°C-40°C 0°C-40°C Operating: Storage:3 -40°C to 70°C -40°C to 70°C

Weight

(with extra battery

option): 12kg (26.4lbs) 12kg (26.4lbs) Size (HxWxD): 145mm x 320mm x 145mm x 320mm x

> 346mm 346mm 5 5/8" x 12 1/2" x 5 5/8" x 12 1/2" x

13 5/8" 13 5/8"

#### NOTES

- All values measured after a one-hour warm-up period.
- Fully floating relative to earth ground. 2
- To prevent damage, batteries should be fully charged before subjected to temperature extremes.

In keeping with our commitment to continuous improvement, ILX Lightwave reserves the right to change specifications without notice and without liability for such changes.

#### ORDERING INFORMATION

LDX-3620 Ultra-Low Noise Current Source

LDX-362038 Extra Battery Option

400164 Small Replacement Battery (1.2A-hour) 400165 Large Replacement Battery (4.5A-hour) CC-305S Current Source/ Laser Diode Mount

Interconnect Cable

CC-306S Current Source/Unterminated

Interconnect Cable

RM-132 Single Rack Mounting Kit

For information call 1-800-459-9459 International Inquiries: 406-556-2481

email: sales@ilxlightwave.com

